

Omni Reaction Stations and Digi-mantle Quick Setup Guide


Set up

1. Place the Dry Block or Mantle into the Omni Station ensuring the connection plug and socket align. **(figure 1)**
2. Place the Omni in the place you would like it and insert the mains leads into the socket at the back of the Omni Station. **(figure 2)**

NOTE: Please ensure that the correct voltage is used.

3. Place glassware into the mantle or into each position on the reaction station.
4. (If required). Fit reflux unit over the dry block with the glassware protruding through the ports. Fit reflux caps and connect the purge tubing.

NOTE: Please refer to the operator guide for reflux use.

5. Turn on the Omni by pressing the I/O switch on the front of the unit. After 'Power-on' the keyboard will illuminate and the front screen will perform a 'self-check' before displaying the module type fitted. **(figure 3)**
6. Press the  key to enter the main menu. The LCD display menu on the front screen will show the current operating condition. **(figure 4)**

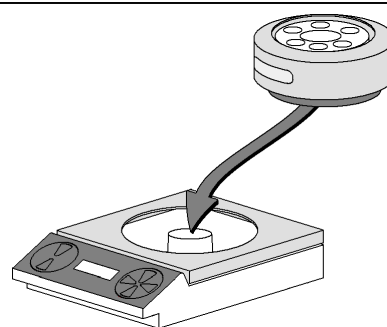


Figure 1

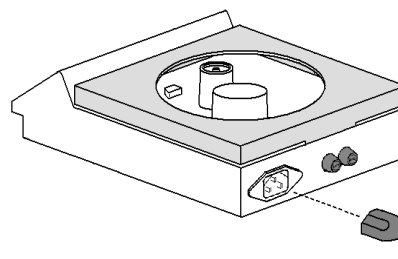


Figure 2

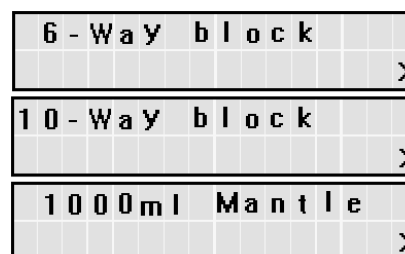


Figure 3

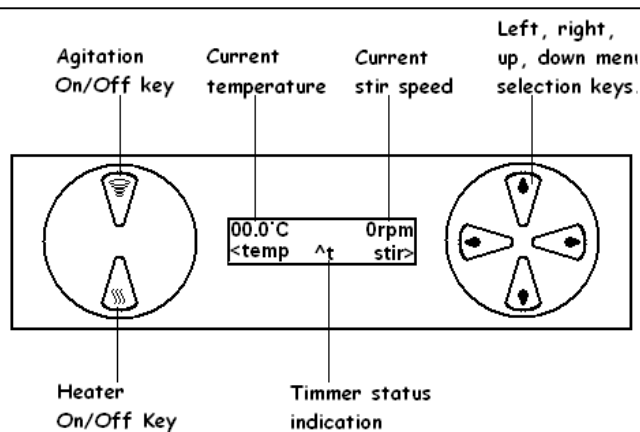


Figure 4